



Improved Productivity with the Niagara Rapid Rinse: A Simple, Low-Cost Solution for Greater Sample Throughput

Authors: Justin Masone and Ryan Brennan

Most ICP and ICP-MS systems incorporate an autosampler. With these systems, there is a significant delay between the time when the autosampler probe enters the sample and the time when the sample reaches the nebulizer. There is a similar delay between the time when the probe enters the rinse and the time when the rinse reaches the nebulizer. If these delays could be eliminated, the analysis time per sample would be reduced. This is exactly what the Niagara CM Rapid Rinse (Niagara) achieves.

In this article, we compare the sample cycle time of a standard ICP-MS system with the sample cycle time of the same ICP-MS coupled with the Niagara, **where a productivity increase of 91% was achieved.**

Introduction

The Niagara is the simplest, lowest-cost method of improving throughput on an existing ICP or ICP-MS system. Increased throughput is achieved by rinsing the nebulizer and spray chamber the instant the sample measurement is completed and continuing to rinse until the next sample has reached the nebulizer. Thus, the rinse is carried out during the time it takes for the sample and the rinse solutions to flow from the autosampler probe to the nebulizer. Typical time saved with the Niagara is around 30%, and, since the valve switches immediately after the read is complete, the amount of sample introduced to the ICP-MS is minimized.¹ This feature drastically reduces cost of consumables and frequency of cleaning, particularly for challenging sample matrices.²

In **Figure 1**, we show a flow diagram of the Niagara in both the Home and Rinse positions. The valve is in the Home position in **Figure 1a**, where the sample is being directed to the nebulizer while the rinse solution is directed to waste. At the completion of the sample measurement, the valve switches instantly to Rinse position and the rinse solution is directed to the nebulizer, as shown in **Figure 1b**, allowing for the nebulizer and spray chamber to begin rinsing immediately, with no uptake delay. The valve remains in the Rinse position while the autosampler probe moves to the next sample. Only when the next sample has made its way through the full length of the uptake tubing does the valve switch back to the Home position (**Figure 1a**).

Figure 1a. Sample to nebulizer; rinse to waste (inject position)



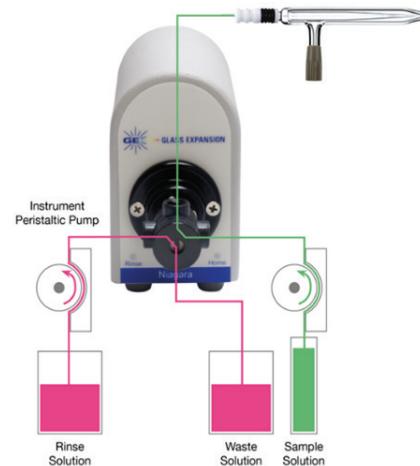


Figure 1b. Rinse to nebulizer; sample to waste (rinse position)

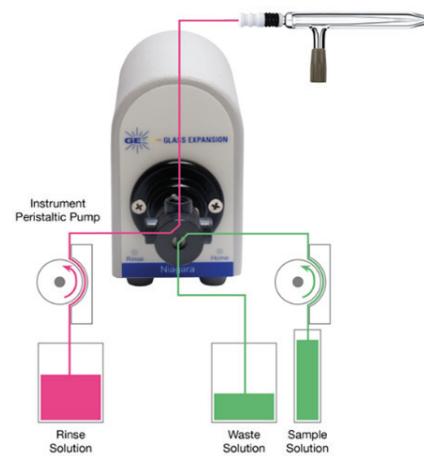

Part of the simplicity in installing and optimizing the Niagara Rapid Rinse lies in the software package. The software includes a Method Wizard (**Figure 2**), which features an Uptake Delay Calculator, Rinse Tubing Calculator, and Time in Sample (TIS) Calculator. The addition of these calculators greatly simplifies method development so that you can quickly add the Niagara to your existing ICP method and optimize for maximum savings.

Figure 2. Niagara Rapid Rinse Method Wizard

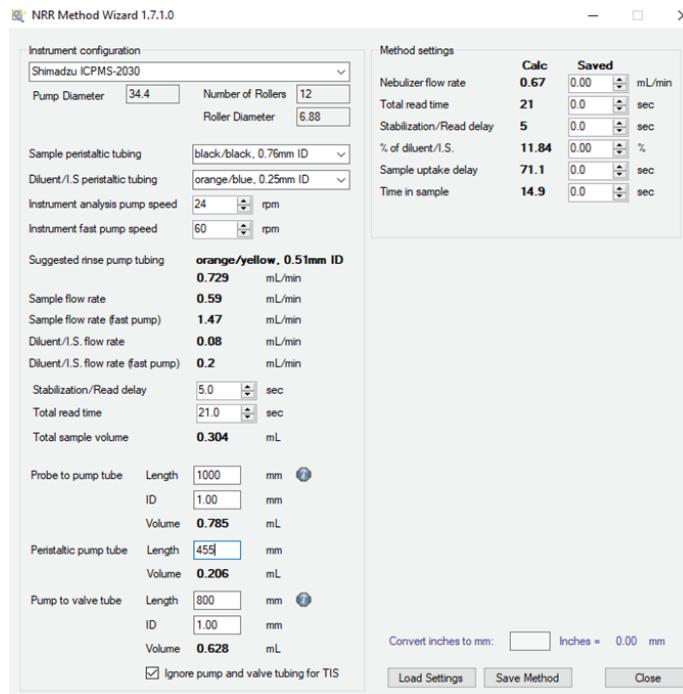

The Uptake Delay Time is the time required for the sample to reach the valve and pass through to waste, ensuring that the first bit of sample diluted with the previous rinse is flushed out to waste. Once the Uptake Delay Time has finished, the Niagara valve will switch from Rinse to the Home position, directing the sample to the nebulizer. The Uptake Delay Time is calculated based on the ID and length of the autosampler probe; peristaltic pump tubing ID and length; and RPM of the peristaltic pump (Fast Pump Speed).

The Rinse Tubing Calculator calculates the optimum peristaltic pump tubing to be used in order to achieve the same flow rate that is used during analysis while using a fast pump during the uptake and post-acquisition rinsing steps of the analysis.

Time in Sample (TIS) is the time required for the autosampler probe to remain in the sample tube during the acquisition. The TIS is based on the total read time, RPM of the peristaltic pump (Analysis Pump Speed) and dimensions of the uptake tubing referenced above.

To improve washout, it is well known that adding a bubble or series of bubbles can help “scrub” the uptake line.¹ However, air bubbles introduced into the plasma can lead to instability and require longer stabilization delay times. With the Niagara, we can easily add a series of air bubbles with our post wash command, without the worry of introducing any air or extra argon into the plasma. To achieve this, the Niagara software sends the autosampler probe to the rinse station after the TIS has completed, followed by a series of commands to slowly raise and lower the probe in the rinse station.

This creates a series of air gaps between the rinse solution, helping to rinse the probe, uptake tubing and sample peristaltic pump tubing. Once the ICP method rinse begins, the air bubbles and any excess sample are directed to waste, while rinse solution is directed to the nebulizer.

Results

A Shimadzu® ICPMS-2030 with a Shimadzu® AS-10 autosampler was used for this study. The standard sample introduction system includes a MicroMist™ nebulizer (0.6 mL/min), a Twister™ spray chamber, and a quartz mini torch, all supplied by Glass Expansion. The Niagara Rapid Rinse accessory was installed using a custom bracket, in order to position the valve as close to the nebulizer as possible, as seen in **Figure 3**.

Figure 3. Glass Expansion's Niagara Rapid Rinse accessory installed on a Shimadzu® ICPMS-2030



Calibration and check standards were created at concentrations typically used in the cannabis industry, and a method was generated for the analysis of the big four: As, Cd, Hg, and Pb. These values are given in **Table 1**. Internal standardization was accomplished for all standards, samples, and blanks automatically using the Glass Expansion Trident™ in-line addition kit. The analysis conditions were optimized for speed and accuracy, and these parameters are shown in **Table 2**. The analysis conditions were not changed when the Niagara was added, as the Niagara has no effect on the analysis portion of the sequence.

Table 1. Calibration standards and analyte registration

Analyte	Internal Standard	ug/L	ug/L	ug/L
⁷⁵ As	¹¹⁵ In	0.0 ppb	1.0 ppb	10.0 ppb
¹¹¹ Cd	¹¹⁵ In	0.0 ppb	1.0 ppb	10.0 ppb
²⁰² Hg	²⁰⁹ Bi	0.0 ppb	1.0 ppb	10.0 ppb
²⁰⁸ Pb	²⁰⁹ Bi	0.0 ppb	1.0 ppb	10.0 ppb

Table 2. Operating conditions.

Parameter	Value
Plasma Stand	
RF Power	1.2 kW
Sampling Depth	5.0 mm
Plasma Gas	8.0 L/min
Aux. Gas	1.1 L/min
Carrier Gas	0.7 L/min
Spray Chamber Temp.	5°C
Collision Cell Gas (He)	6.0 mL/min
Analysis Conditions	
Number of Scans	20
Number of Replicates	3
Integration Time	1.0 s

The standards and samples were first analyzed without the Niagara installed, where the sample-to-sample time was 160 seconds, which translates to 22 samples per hour.


Following the default configuration, the Niagara was installed and its performance optimized utilizing the Method Wizard as referenced above. With the Niagara installed, the rinse time, as programmed in the Shimadzu® LabSolutions software, was set to 0, as opposed to 75 seconds in the default configuration. The sample-to-sample time with the Niagara installed was 85 seconds, which translates to 42 samples per hour. This results in a productivity improvement of 91%, allowing an extra 160 samples analyzed per 8-hour day, or an extra 41,760 samples analyzed per year (8h/d, 5d/w).

Calibration curves for the big four analytes, with the Niagara installed, are presented in **Figure 4**, which highlight the accuracy of the Niagara system.

Element/Mass information

Element	Mass	Type	Cond	Int time	Int Mode	IntStd
As	75	QUANT	DBG	1.0	Auto	In 115 (DBG)
Bi	209	INT	DBG	1.0	Auto	
Cd	111	QUANT	DBG	1.0	Auto	In 115 (DBG)
Hg	202	QUANT	DBG	1.0	Auto	Bi 209 (DBG)
In	115	INT	DBG	1.0	Auto	
Pb	208	QUANT	DBG	1.0	Auto	Bi 209 (DBG)

Figure 4a. Calibration curve for As

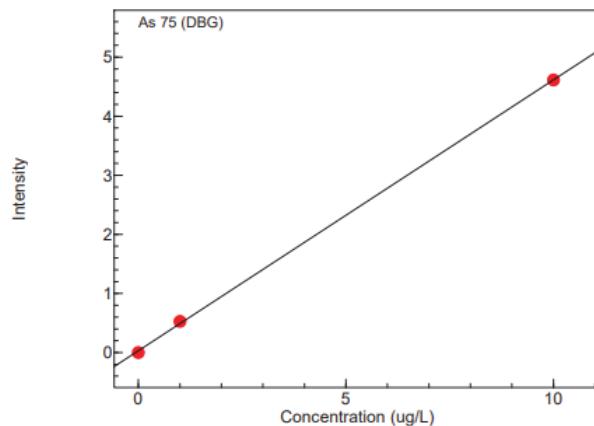


$$\text{Conc.} = 2.181071 * I - 0.0706261$$

r = 0.99991 BEC = 0.0706261 (ug/L)
Weight: None Origin: None


Class	Sample Name	Set Conc.	Intensity	Conc.	Diff of Conc.	Excl
CAL1	blk	0.0000000	0.0007319	-0.0690	-0.0690	
CAL2	1 ppb	1.0000000	0.5265939	1.08	0.0779	
CAL4	10 ppb	10.00000	4.613213	9.99	-0.0089	

Figure 4b. Calibration curve for Cd

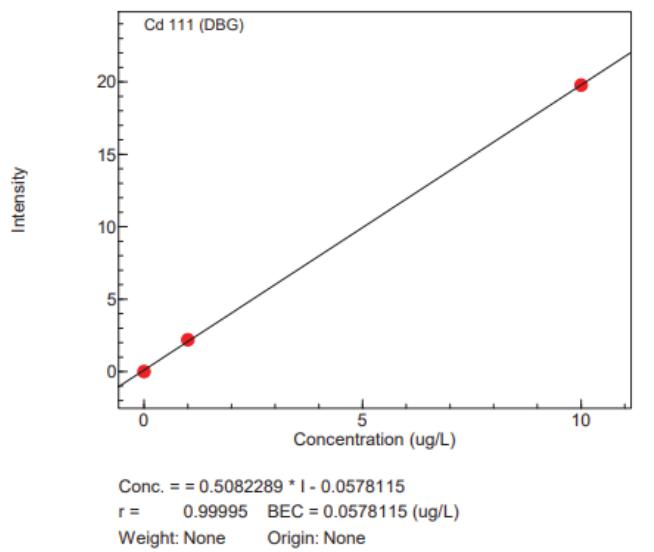


Class	Sample Name	Set Conc.	Intensity	Conc.	Diff of Conc.	Excl
CAL1	blk	0.0000000	0.0076362	-0.0539	-0.0539	
CAL2	1 ppb	1.0000000	2.200724	1.06	0.0607	
CAL4	10 ppb	10.000000	19.77669	9.99	-0.0067	

Figure 4c. Calibration curve for Hg

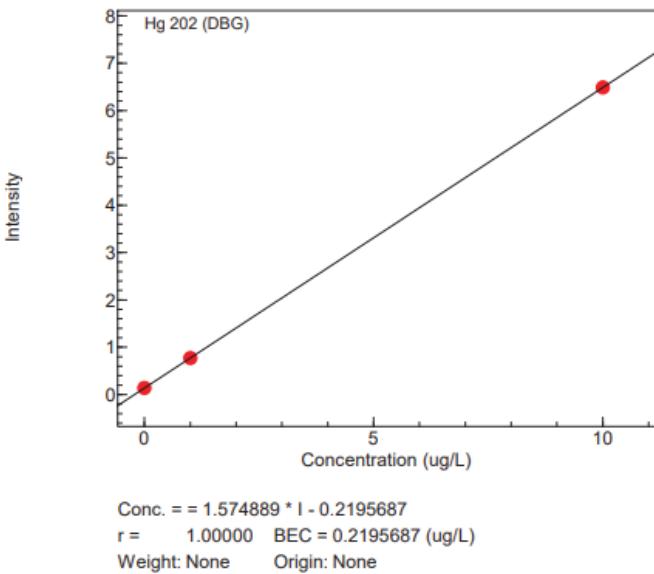

Class	Sample Name	Set Conc.	Intensity	Conc.	Diff of Conc.	Excl
CAL1	blk	0.0000000	0.1412768	0.0029	0.0029	
CAL2	1 ppb	1.0000000	0.7723206	0.997	-0.0032	
CAL4	10 ppb	10.000000	6.489277	10.0	0.0003	

Figure 4d. Calibration curve for Pb

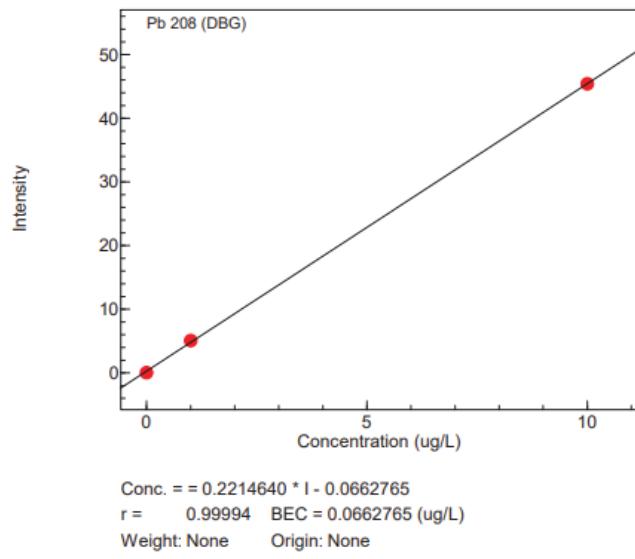

Class	Sample Name	Set Conc.	Intensity	Conc.	Diff of Conc.	Excl
CAL1	blk	0.0000000	0.0510705	-0.0550	-0.0550	
CAL2	1 ppb	1.0000000	5.093903	1.06	0.0618	

Conclusion

The Niagara is the simplest, lowest-cost method of improving throughput on an existing ICP or ICP-MS system. The results from the study show that, once the Niagara Rapid Rinse is installed and optimized, significant time savings are achieved. This is done by rinsing the nebulizer and spray chamber the instant the sample measurement is completed and continuing to rinse until the next sample has reached the nebulizer. Thus, the rinse is carried out during the time it takes for the sample and the rinse solutions to flow from the autosampler probe to the nebulizer. While typical time savings are around 30%, in this particular case, time savings of 47% were achieved, which would result in an extra 41,760 samples per year.

References

1. Glass Expansion June 2016 Newsletter, *A Simple Low Cost Approach to Improving ICP Productivity*.
2. Glass Expansion June 2017 Newsletter, *Considerations when Analyzing Real-World Samples by ICP Spectrometry*.